

HITACHI
Inspire the Next



STF3
Tube Sealer

himac
centrifuges

Tube Sealer model STF3

The STF3 is heat-welding tube sealer to use the seal tubes with fixed angle rotors, vertical rotors and neo angle rotors.

Features

- Simple & Easy Operation
 1. Set seal tubes in the appropriate tube rack then set the tube rack on the rack guide of the STF3.
 2. Adjust position of the tube rack in order the tube inlet comes under a heater of the STF3.
 3. Pull the handle down to contact the heater the tube inlet and hold the handle for 1 to 2 seconds. Then pull the handle down to the end and wait until the HEAT Lamp goes off. (It takes about 40 seconds.)
- CE Marking is qualified.
- HEAT Lamp indicates the status of the heater.
- Tube racks are compatible with our former model STF2.



Specifications

Model	STF3 (110-120V equipment)	STF3 (200-240V equipment)
Seal Method	Direct welding by the molding heater	
Standard	CE marking, cCSAus	CE marking
1. EMC	EN61326-1	
2. Product Safety	EN61010-1	
Dimension (mm)	132 (W) x 225 (D) x 260 (H) < when raising the handle : 320 (H) >	
Weight (kg)	8.0	
Power	99 - 132 VAC, 2A, 50/60Hz	180 - 264 VAC, 2A, 50/60Hz
Part No.	90132501	90132502

Tube Rack

Please order appropriate tube rack*1 that is suitable for the seal tube to be used.

Part No.	Description	Applicable Tube	Remarks
S201778G	Tube Rack (G2)	1.5PA seal tube	
S201778F	Tube Rack (G)	2PA seal tube	
S201778E	Tube Rack (B2)	3.5PA seal tube	
S201778H	Tube Rack (B3)	4PA seal tube	
S201778A	Tube Rack (B)	5PA seal tube	
S201778J	Tube Rack (B4)	6.5PA seal tube	
S201778L	Tube Rack (C2)	8PA seal tube	
S201778B	Tube Rack (C)	12PA seal tube	
S201778C	Tube Rack (E)	40PA seal tube	
S201778M	Tube Rack (F2)	3.5PA seal tube	Not for STF1
S201778K	Multi Rack	2, 4, 5, 6.5, 12 and 40PA seal tube	Not for 1.5, 3.5, 8, and 94PA seal tube

*1 : sold separately

The photos used on this leaflet are STF3 (200-240V equipment).



CAUTION:

For safe and proper use of your machine, carefully read and follow the instructions in the instruction manual.

- All specifications are subject to change without advanced notice.
- Actual color may vary from the color of the photos on this catalogue, due to printing condition.
- Due to safety reason, installation environments, operating environments and conditions may be restricted.
- Unless specially mentioned, products and/or operation panel of the photos are standard specification.
- For further information, please contact your nearest Hitachi Koki's representative.

Manufactured by :

Hitachi Koki Co., Ltd.

Life-Science Instruments Division

1-8-12 Shiba-kouen, Minato

Tokyo 105-0011 JAPAN

TEL : +81-3-5733-0206

FAX : +81-3-5733-0207

URL : <http://centrifuges.hitachi-koki.com/>